Smartphone Microscope

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Dear customer,

Thank you for purchasing our product. Please read the following instructions carefully before first use and keep this user manual for future reference. Pay particular attention to the safety instructions. If you have any questions or comments about the device, please contact the customer line.

www.alza.co.uk/kontakt

+44 (0)203 514 4411

Importer

Alza.cz a.s., Jankovcova 1522/53, Holešovice, 170 00 Prague 7, www.alza.cz

Thanks for choosing our Smartphone Microscope. Please read and keep instructions carefully to ensure proper functions of your new Microscope.

Packaging List

1. Smartphone Microscope
2. Multi-function Stretch Clip
3. Type-C Charging Cable
4. Cleaning cloth
5. EVA Pads (2/3/4mm)
6. User Manual




Microscope Intro


Operation

Multi-functional Stretch Clip Installation

Gently push your phone into the stretch clip, and make sure the 17mm-thread port is center align with main lens of the phone.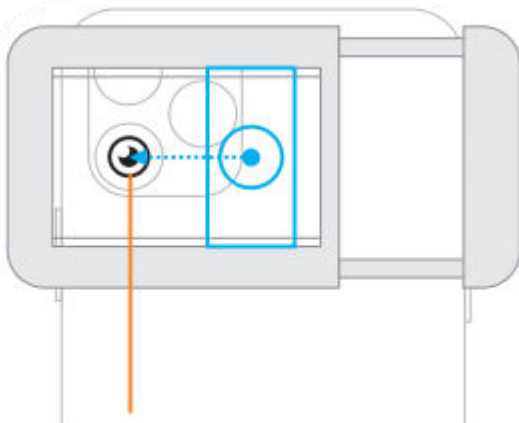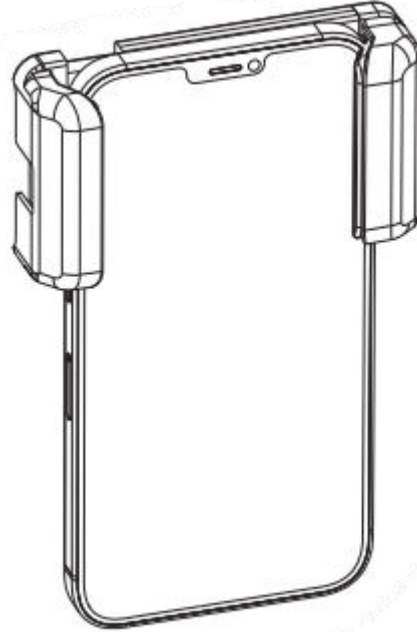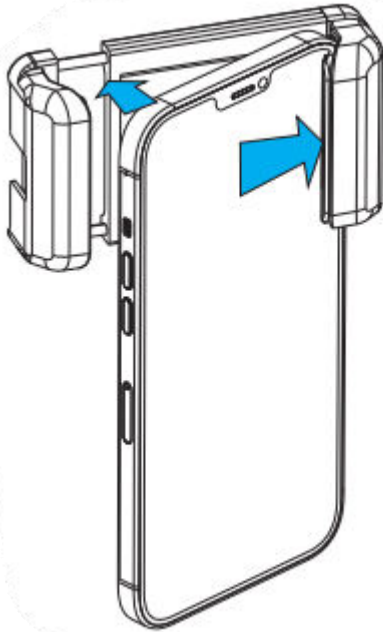


Then slide the clip's metal plate and align the 17mm-thread port with the center of smartphone's main camera.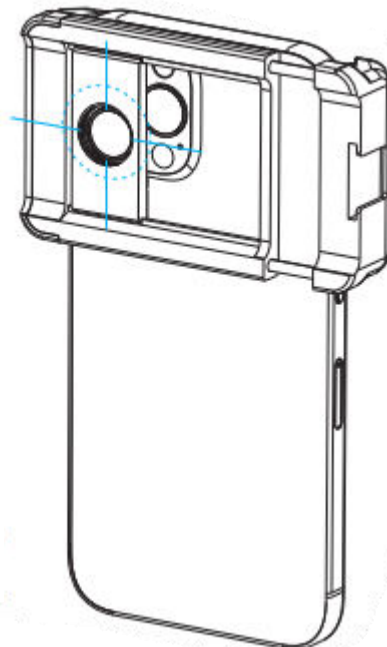

Operation


Microscope Installation

Screw the microscope clockwise into 17mm threaded port of the clip. Then press the power switch to turn on LED's.
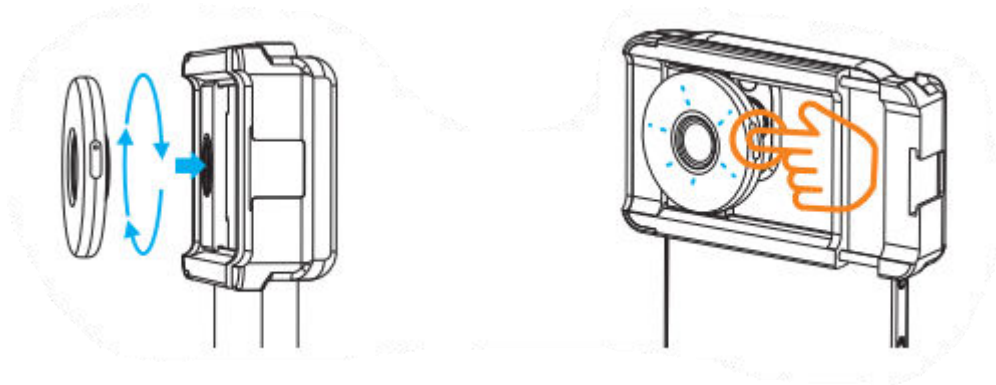


Microscope with Smartphone Shooting

Please keep the same smartphone and the microscope in horizontal or vertical direction for shooting. The best focusing distance is 5.4mm to 6.8mm.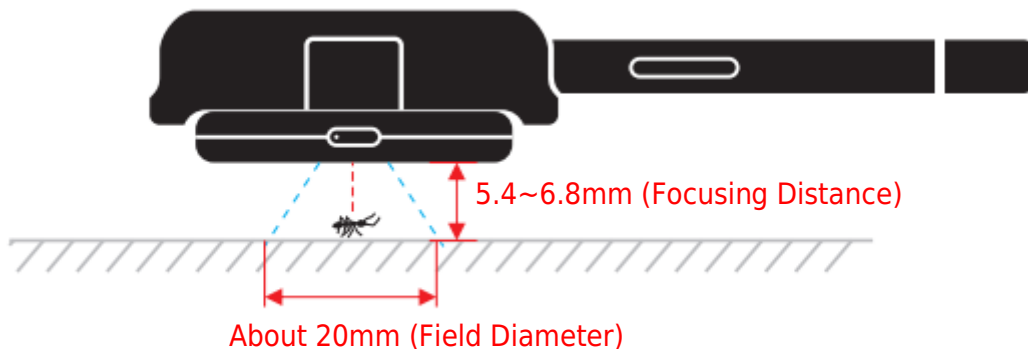


- Please get focused on the subject before shooting.
- Please do not tilt the phone and microscope in during shooting.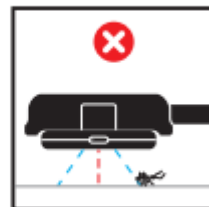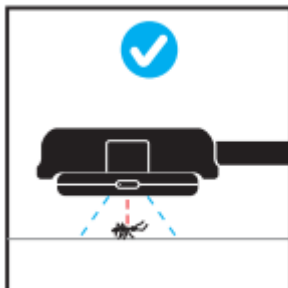

Microscope Charging Instructions

Connect cable and type-c charging Port, and plug other side into the power (Computer/Power Bank/USB charger).


- The red light is on when it's charging, then the green light is on when the charge is completed.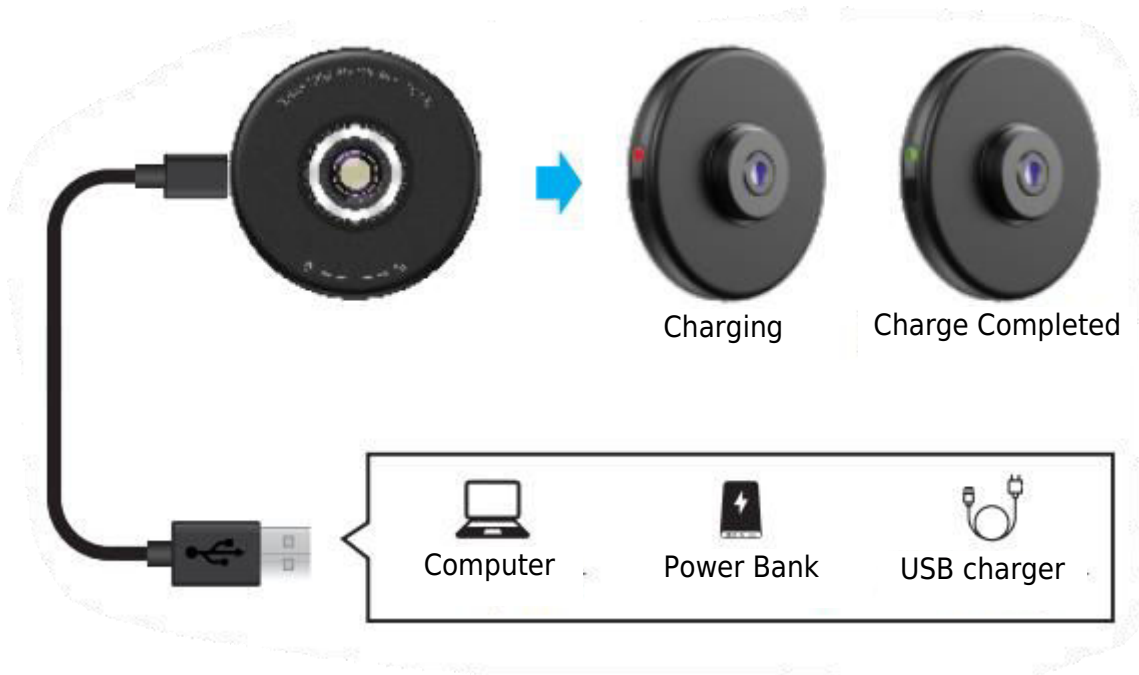

Note:


- The microscope charging time is about 50 minutes, and using time is about 120~150 minutes.
- Please do not charge in microscope shooting time.
- To keep the battery alive, please charge the microscope every three months if you don't use it for a long time.
- When charging, do not let the interface be invaded by water or small debris, otherwise the interface may easily cause short circuit, leakage and other faults.

Multifunctional Stretch Clip Introduction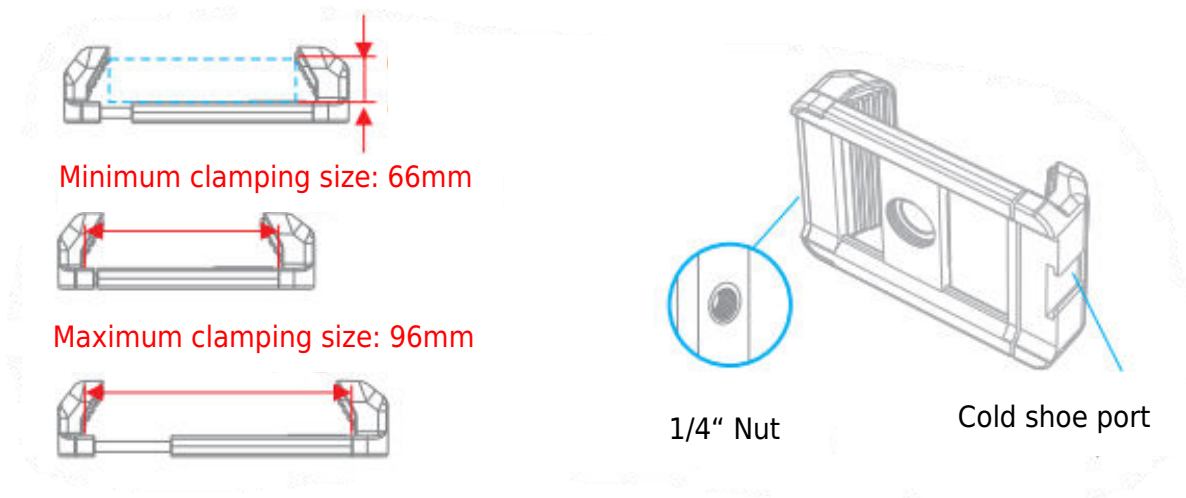


Pay attention to the installation direction of the clip

If the clip installation direction is not suitable, the 17mm thread port will not be aligned with the phone's main camera, please reverse the direction of the clip and then install it.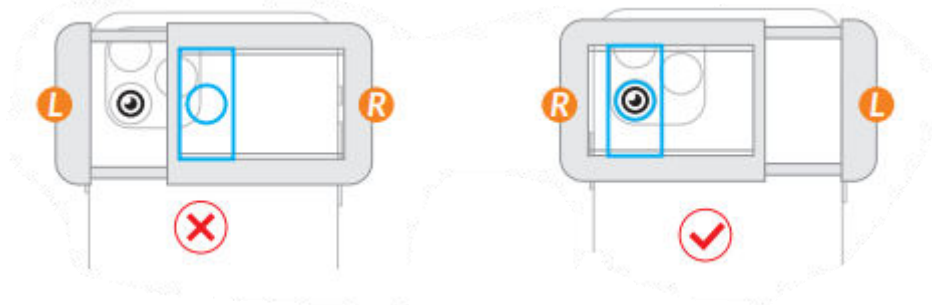

Included EVA Pad

If your smartphone camera is raised, please choose EVA pads to stick on clip inner (the pads are available in 3 thickness: 2mm, 3mm, 4mm) to fill the gap on your phone back to make it flushes with the camera.

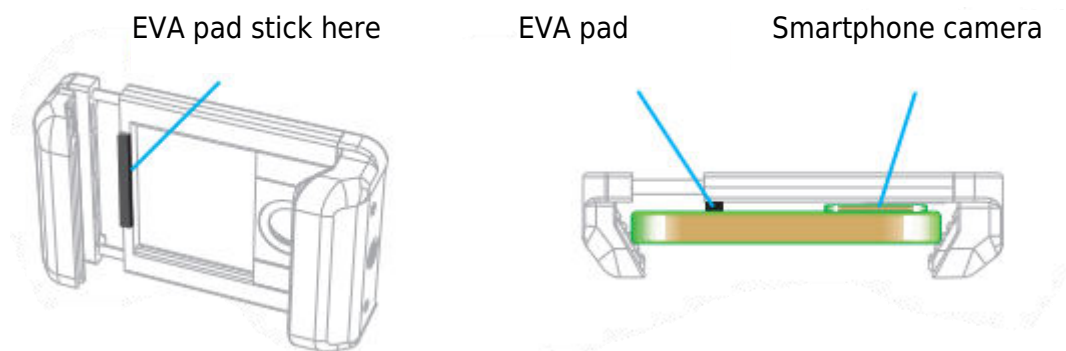Warranty Conditions

A new product purchased in the Alza.cz sales network is guaranteed for 2 years. If you need repair or other services during the warranty period, contact the product seller directly, you must provide the original proof of purchase with the date of purchase.

The following are considered to be a conflict with the warranty conditions, for which the claimed claim may not be recognized:

- Using the product for any purpose other than that for which the product is intended or failing to follow the instructions for maintenance, operation, and service of the product.
- Damage to the product by a natural disaster, the intervention of an unauthorized person or mechanically through the fault of the buyer (e.g., during transport, cleaning by inappropriate means, etc.).
- Natural wear and aging of consumables or components during use (such as batteries, etc.).
- Exposure to adverse external influences, such as sunlight and other radiation or electromagnetic fields, fluid intrusion, object intrusion, mains overvoltage, electrostatic discharge voltage (including lightning), faulty supply or input voltage and inappropriate polarity of this voltage, chemical processes such as used power supplies, etc.
- If anyone has made modifications, modifications, alterations to the design or adaptation to change or extend the functions of the product compared to the purchased design or use of non-original components.

EU Declaration of Conformity

Identification data of the manufacturer's / importer's authorized representative:

Importer: Alza.cz a.s.

Registered office: Jankovcova 1522/53, Holešovice, 170 00 Prague 7

CIN: 27082440

Subject of the declaration:

Title: Smartphone Microscope

Model / Type: MS009

The above product has been tested in accordance with the standard(s) used to demonstrate compliance with the essential requirements laid down in the Directive(s):

Directive No. 2014/30/EU

Directive No. 2011/65/EU as amended 2015/863/EU

WEEE

This product must not be disposed of as normal household waste in accordance with the EU Directive on Waste Electrical and Electronic Equipment (WEEE - 2012/19 / EU). Instead, it shall be returned to the place of purchase or handed over to a public collection point for the recyclable waste. By ensuring this product is disposed of correctly, you will help prevent potential negative consequences for the environment and human health, which could otherwise be caused by inappropriate waste handling of this product. Contact your local authority or the nearest collection point for further details. Improper disposal of this type of waste may result in fines in accordance with national regulations.

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